


Installation of NucleoView 2.x on Windows 7 operating system (32 bit only)


This document will also apply for installation of the USB Driver for SemenView and SomaticView applications.

In order to perform an installation of the USB driver on Windows 7 (32 bit version), the following message will most likely appear.

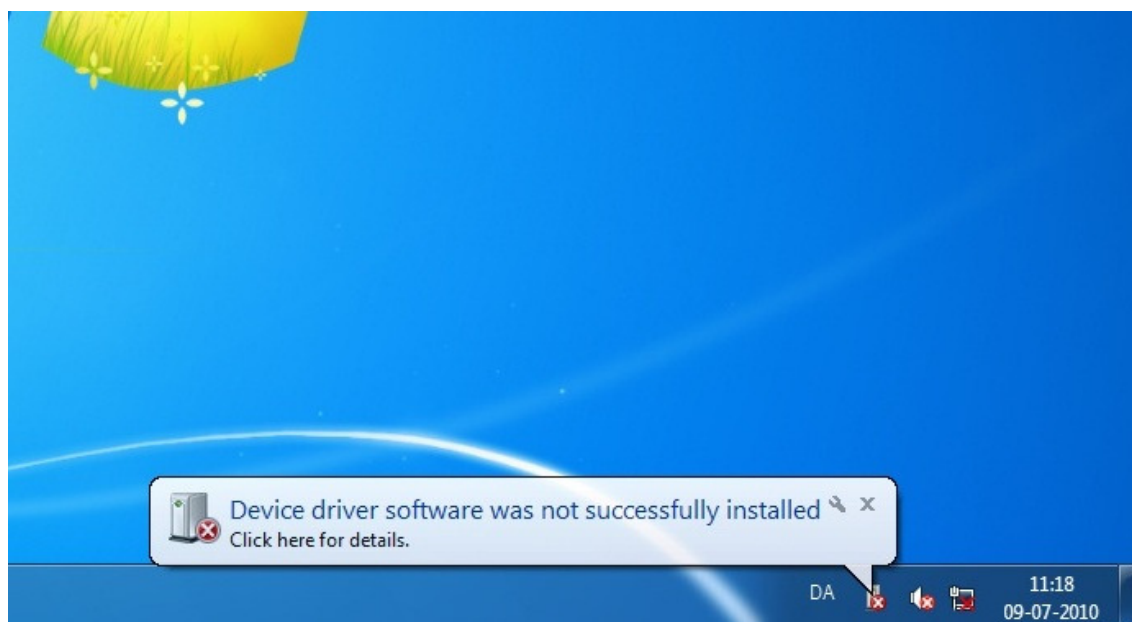

This message tells that Windows 7 not automatically could install the USB driver for the NucleoCounter Instrument. Please ignore this message and follow the instructions hereafter.

Due to different user levels and security settings of Windows 7, we recommend that the local IT administrator supports in this matter.

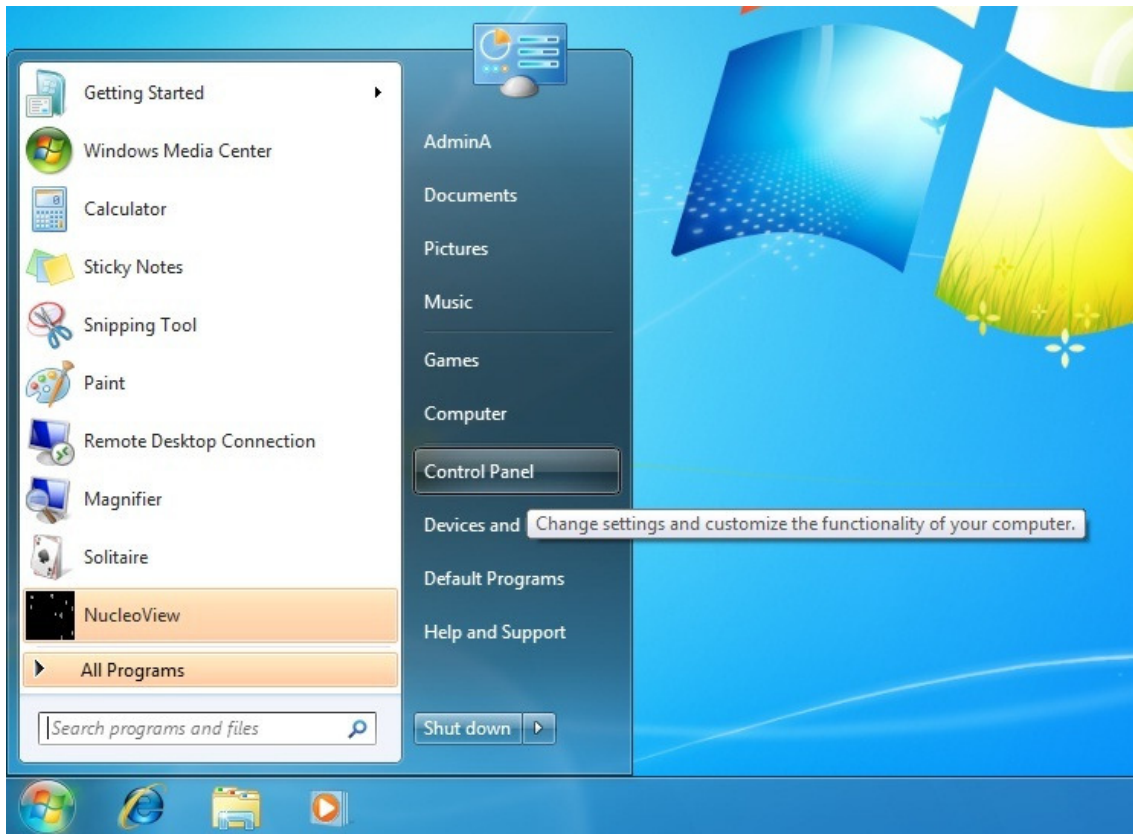


Through the Start button (lower left corner typically) activate the “Control Panel”



Select “System and Security”




At the "System" menu select "Device Manager"

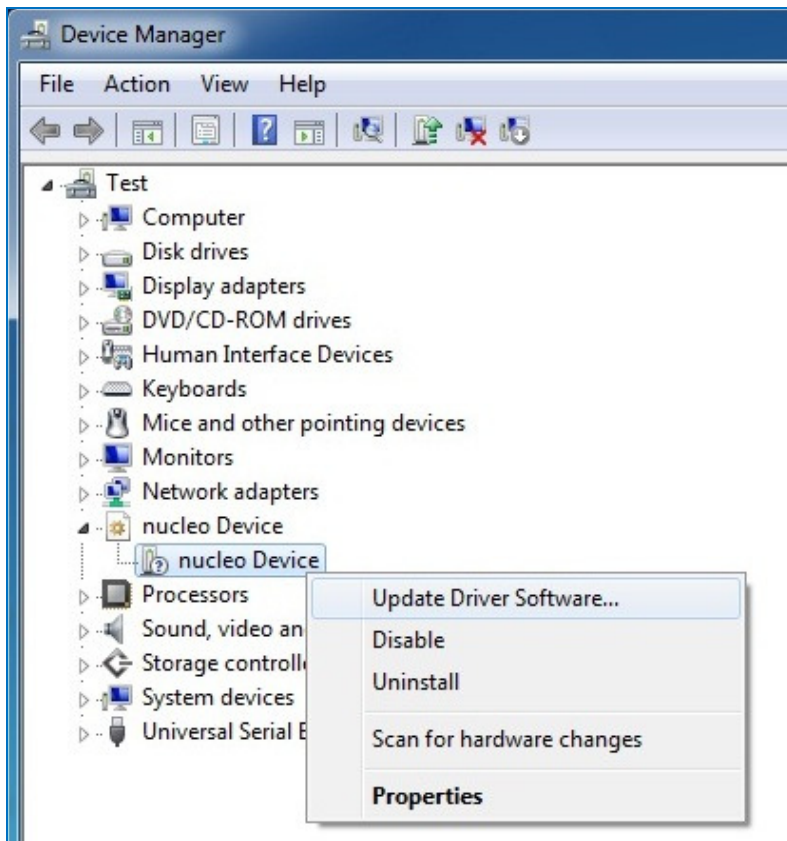

Locate and open the "nucleo Device" item. Click on the right mouse button and select "Update Driver Software"



Select "Browse my computer for driver software"

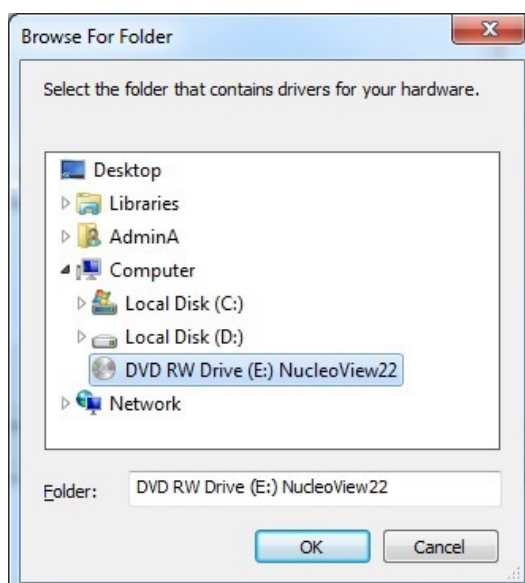

In the "Browse for Folder" dialog locate the drive with the "NucleoView22" CD-ROM¹, click OK and in a moment the system will report a successful driver installation. If not, contact the local IT administrator and check for the following: Is a 32 bit operating system present? Are appropriate rights present for installation of new drivers?

Otherwise please contact ChemoMetec on +45 48 13 10 20 and ask for assistance, or send an e-mail to contact@chemometec.com.

¹ Or SemenView or SomaticView CD-ROMs